



Manual update of Octoview software

Document-Version 1.1 09/2025

Author: Dr. rer. nat. Patrick Vogel

eMail: Patrick.Vogel@phase-vision.com

All rights reserved. Copyright 2025 by phase VISION GmbH





1. Manual Octoview update

Step 1: Unpack the provided zip-file consisting of the new files:

File structure

|-- **OctoView**

 |-- **Win64**

 |-- **Release**

 OctoView.exe

 MPI_server.exe

 Updater.exe

 libopenblas.dll

 ChangeLog.html

 pd_icon.png

 _release.docx

 OctoView_features.docx

 |-- **Documentation**

 Documentation.pdf

Step 2: Go to the installation folder

default:, C:\Program Files\Pure Devices\PD OctoView\

Step 3: Replace all files with the new ones.

HINT: Before replacing the old files, they can be renamed for backup purposes.

Step 4: Finalization

Check **_release.docx** for new features and updates

Check **documentation.pdf** for new updates.